



BPE

Bovine Pituitary Extract

Catalog Number: S-011-1

1 ml

Instructions for storage and use

Product Description

Bovine pituitary extract (BPE) is a sterile solution containing approximately 13 mg/ml total protein.

Intended Use

BPE is intended for use in conjunction with an appropriate basal medium and other additives for the routine culture of human epithelial cells. ***This product is for research use only. Not for use in animals, humans, or diagnostic procedures.***

Caution: If handled improperly, some components of this product may present a health hazard. Take appropriate precautions when handling this product, including the wearing of protective clothing and eyewear. Dispose of properly.

Storage and Stability

BPE is stored at -20° C at our facility and is shipped on dry ice. Upon receipt, the product should be stored at -20° C in a freezer that is not self-defrosting. When stored at -20° C the product is stable until the expiration date shown on the label.

After long-term storage at -20° C, BPE may contain a small amount of precipitate. This precipitate is formed from cold-insoluble material in BPE and will not affect the performance of the product.

Thawing

To thaw, place the product in a 37° C water bath or overnight at 4° C. If thawed in a water bath, do not leave the product at 37° C after the product has thawed.

Terms and Conditions of Sale

Cascade Biologics, Inc. (hereinafter, CBI) warrants that its products will perform according to the information provided in various publications that it distributes, and as described herein for the intended shelf life of the product when stored under the conditions prescribed by CBI. If any product does not perform according to the published information provided by us, CBI will replace the product free of charge to the original customer. The remedy of product replacement shall be the customers' sole and exclusive remedy for defective product, unless CBI is unable to deliver replacement product, in which case CBI shall reimburse the customer for the purchase price of the defective product. Customer understands that the foregoing limited warranty is in lieu of all other warranties and CBI hereby disclaims all other warranties including, but not limited to, implied warranty of merchantability or fitness for adequacy for any particular purpose or use. CBI shall not be liable to customer or to any party claiming through customer for any incidental or consequential damages, including, but not limited to any lost profits, lost savings, or lost business, whether arising out of contract, tort, or otherwise. Customer acknowledges that products are intended for research use only and are not to be used in animals, humans, or diagnostic procedures.